

# Onto the trinocular tube

For all Motic trinocular microscopes we have several c-mount adapters available. The c-mount adapter needs to be chosen according to the chip size of the camera. Each Moticam (excluding the gooseneck cameras) comes standard with a CS/C-mount connection; the c-ring is included in the package.

Just fix the c-mount adapter on the trinocular port and connect the Moticam to the c-mount adapter. This combination will give you the freedom to continue observing with both eyes, while having the Moticam connected.

Our Moticam cameras can also be used on non-Motic microscopes, via the universal c-mount connection. Just remember that you will need in this case the c-mount adapter from the manufacturer of your microscope.

